

Operating Instructions



863 Valley View Road, Eighty Four PA 15330 USA
Tel: 724-941-9701 Fax: 724-941-1369 e-mail: skctech@skcinc.com

Tedlar® Sample Bags Catalog No. 231-01-TCLP

The 231-01-TCLP is a disposable 1-liter, Tedlar film sample bag with a single stainless steel fitting designed for use with a Zero Headspace Extractor for the EPA TCLP method. The use of Tedlar sample bags with stainless steel fittings has been validated by the EPA.

Operation

The bag is designed to be attached directly to the extractor vessel via a special stainless steel coupler (SKC Catalog No. 231-01-ZHE).

To open the fitting, remove the knurled screw cap on the fitting. Attach the coupler to the bag by screwing onto the bag fitting, then attach directly to the ZHE unit.

Note: SKC 231-01-TCLP bags are designed for single use applications and are not intended for the storage and transport of hazardous materials.

© Tedlar is a registered trade name of DuPont Corporation.

Notice: This operating instruction may not address all safety concerns (if any) associated with this product and its use. The user is responsible for determining and following the appropriate safety and health practices and regulatory limitations (if any) before using the product. The information contained in this document should not be construed as legal advice, opinion, or as a final authority on legal or regulatory procedures.

www.skcinc.com

Form # 3783 Rev 1001