



Exciting News!

On July 15, 2019, SKC Inc. acquired Colormetric Laboratories, Inc. (CLI).

SKC is excited to add CLI surface and dermal assessment and control tools to our highly respected, long-standing sampling solutions that help you maintain the health and safety of workers. SKC looks forward to continuing to provide the same great CLI products with excellent customer and technical support. Learn more about SKC Inc. at <https://www.skcinc.com/about>.

Surface SWYPE for Aliphatic Amines Cat. No. 769-1025

1. Lightly spray the area or item (e.g., workbench, tool, control knob) with Developing Solution for Aliphatic Amines Cat. No. 769-1065.
2. Wait at least 30 seconds and then wipe with a Surface SWYPE pad.
3. Allow 3 minutes for the color to develop on the pad. A fuchsia color indicates that aliphatic amines may be present. This color reaction relies on detecting the alkaline pH of aliphatic amines.

SKC Limited Warranty and Return Policy

SKC products are subject to the SKC Limited Warranty and Return Policy, which provides SKC's sole liability and the buyer's exclusive remedy. To view the complete SKC Limited Warranty and Return Policy, go to <http://www.skcinc.com/warranty>.